



How To Change Your Innocage®



IMPORTANT WARNING

It is not recommended to use Innocage[®] IVC Mouse and Innocage[®] IVC Rat outside of the Innorack[®].

Do not leave animals in an unventilated IVC cage for more than 12 hours.



Remove the water bottle from the dirty Innocage[®].

Open the cage by lifting the right front corner.



2



3

Transfer the animals to a clean cage. Be sure to check clean cage is free of dents. If cage bottom is dented, do not use.

Replace the food tray;
refill it with food using the MAX
line on the tray as a
guide.
Do not overfill.





Place the lid on the Innocage®. Secure it by gently pressing down on the corners. Then slide the cardholder into place on the front of the Innocage®.

If using a pre-filled Aquavive® bottle, remove the seal from the cap. Place water bottle in the lid and check for water flow.





7

Return the Innocage[®] in the Innorack[®] - make sure the water bottle is facing the front.

TIP: When cage is docked correctly, you can feel the cones on the top of the cage lock into place in the rack.

CAGE CHANGE IS COMPLETE!



*Did you know
all our cages for mice
and rats are available
pre-bedded with corn
cob or ALPHA-dri®?*

Disposable Cages for Mice:

- Innocage® IVC
- Innocage® IVC Dual Filter
- Innocage® IVC Large Filter
- Innocage® Containment
- Innocage® Static
- Innocage® Transport

Disposable Cages for Rats:

- Innocage® IVC
- Innocage® Static Rat
- Innocage® Tall Rat

For a complete list of Innovive caging and accessories, visit www.innovive.com



WE RECYCLE!

Call Innovive for more details:

(866) 432-2437 or email recycling@innovive.com

Innovive

www.innovive.com

Less Washing, More Science